


Pill Counting

DRX-4C2 Quick Reference Guide

To enter the counting mode, simply press the F button and follow the steps below.



1. Place a container on the pan and tare it by pressing the "T" button.

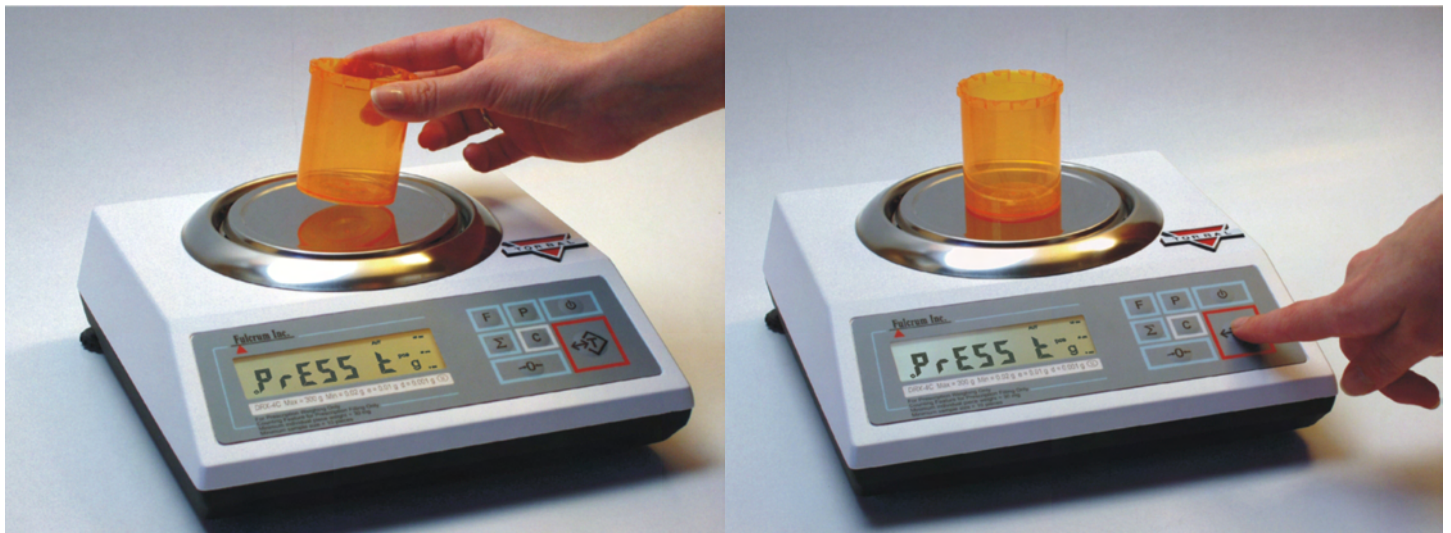

2. Confirm the sample size by pressing the T button.



3. Manually count 10 pills and place them in the container located on the pan.



4. Press the T button to confirm and accept the sample weight.



5. Begin to fill the remainder of the pieces.



For a detailed description of this feature please turn to Chapter 12 of the instruction manual.

For technical questions or assistance please call toll free: 1-866-473-6900

Fulcrum Inc.
 23 Carol Street.
 Clifton, NJ 07014
 (973) 473-6900
 (973) 777-8302

<http://www.fulcruminc.net>